Túi ổn định

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Burstable Pouch Sterile Session Wipes

A complete range of sterile sporicidal, biocidal and detergent wipes in a session wipe.

Pouch compartments maintain wipe and product efficacy for full shelf life.

USE INSTRUCTIONS:

Remove the pouch from the outer dust bag. With the inner bag intact, roll the foil pouch/liquid compartment and apply downward pressure to burst the liquid through the frangible seal. Leave for one minute to saturate the wipes. Remove the foil pouch from the outer bag. Tear open the end of the bag and peel open the foil pouch to gain access to the wipes. Use a recommended wiping technique to clean and disinfect the surface.

TECHNICAL WIPE DATA:

Lint Grade 1. 68gsm Polyester (45%) / Cellulose (55%) material.
230mm x 230mm (9" x 9").
Thickness 0.32mm.

STORAGE:

Store in original packaging away from sunlight and extremes of temperature. Full information on the storage and handling is available on the Material safety Datasheets (MSDS)

SAFE HANDLING:

Always wear gloves and goggles or face protection. Always read the label and MSDS before use.

MATERIAL COMPATIBILITY:

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

BENEFITS:

- HA (Hypochlorous Acid) suitable for the aseptic transfer process.
- Fully validated formulations
- <60 seconds activation time
- In-use shelf life of 10 hours
- Session wipe format minimises waste
- Any combination of AN, QT, OX and HA are ideal rotational partners
- Manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.
- Sterile wipes until the point of use.
- Multiple bags for cleanroom transfer
- Manufactured to GMP
- The solution is filtered through a 0.2 micron filter and aseptically filled into pre-gamma irradiated packaging (25-45kGy) and tested to give a sterility assurance level SAL of 10~6

SPECIFICATION:

The burstable wipes are available in a choice of five presentations: -

OX Sporicidal preparation of Hydrogen Peroxide 6%/Peracetic Acid 0.4% v/v. Rapid action 2 minute kill time. 2 year shelf life.

HA Sporicidal preparation of Hypochlorous Acid. A safe handling, low odour, low residue. Rapid action 1minute kill time. 1 year shelf life.

QT Biocidal quaternary ammonium compound disinfectant cleaner. A rapid action, non-oxidising biocide. 2 year shelf life.

AN Biocidal diamine disinfectant. A broad spectrum, rapid action, low residue, QUAT free. 2 year shelf life.

Biocides/Detergent are diluted with WFI quality water.

InSpec wipes are manufactured to GMP in an ISO 6 cleanroom. The products are filled and sealed in an ISO class 5 environment.

InSpec™ Burstable Pouches			
Product	Pack Size	Case Size	Product Code
InSpec™ HA Burstable Pouch	15 wipes/pouch	20 pch	HAPW50BP15S
InSpec™ OX Burstable Pouch	15 wipes/pouch	20 pch	OXPW40BP15S
InSpec™ AN Burstable Pouch	15 wipes/pouch	20 pch	ANWF120BP15S
InSpec™ QT Burstable Pouch	15 wipes/pouch	20 pch	QTWFI30BP15S


Vietnam Official Distributor:

Decon Solutions Co., LTD

No 62 Road 64, Thanh My Loi Ward, District 2, HCM City, Vietnam
H: 09 0890 1955 F: +84 28 6287 3181 E: info@deconsolutions.com.vn
T: +84 28 6287 3180 - 38 143 585 W: www.deconsolutions.com.vn

Instructions for Use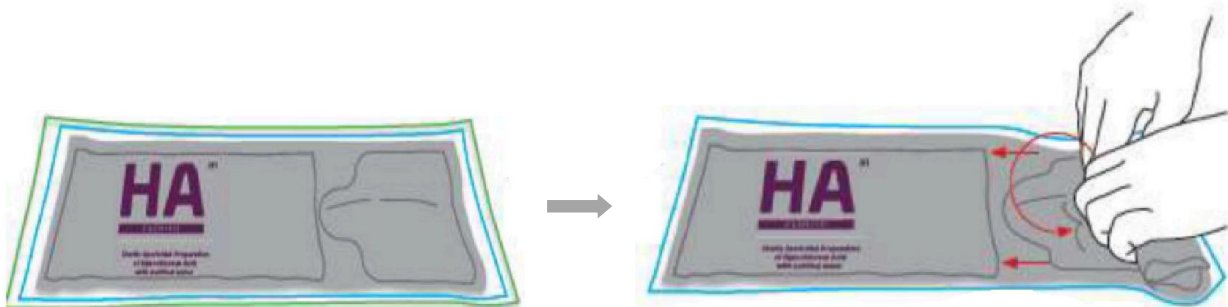

1 With outer packaging still in place roll the pouch from the liquid end and apply pressure to force liquid across the frangible barrier to wet the wipes.

2 With outer packaging still in place roll the pouch from the liquid end and apply pressure to force liquid across the frangible barrier to wet the wipes.


3 Ensure liquid agent is transferred to wipes section.


1MINUTE

4 Wait for minute, then remove the dear outer bag and access wipes.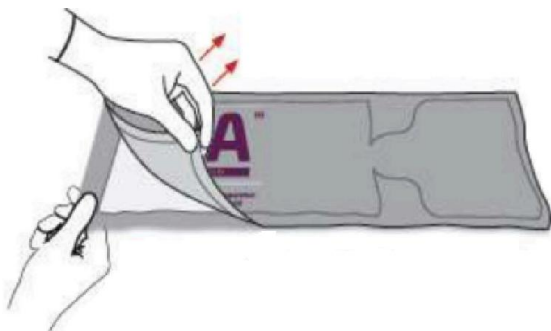

5 To access the wipes, peel the pouch open at the Wipe end of the pouch.

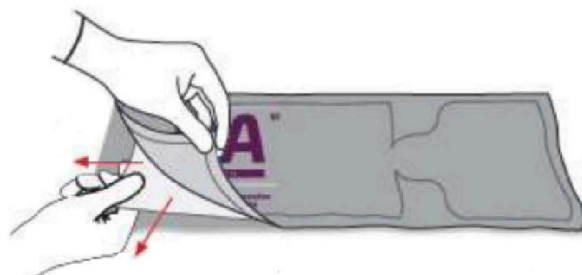6 Use within 10 hrs, re-covering as necessary.

No 62 Road 64, Thanh My Loi Ward, District 2, HCM City, Vietnam
H: 09 0890 1955 F: +84 28 6287 3181 E: info@deconsolutions.com.vn
T: +84 28 6287 3180 - 38 143 585 W: www.deconsolutions.com.vn

Vietnam Offical Distributor:

Decon Solutions Co., LTD