# DE

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# DE

## Sterile Denatured Ethanol 70% v/v with WFI Quality Water

Low Residue, Rapid action, Fast Drying, Alcohol Disinfectant

**USE INSTRUCTIONS:**

InSpec DE is designed for spraying, wiping and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. Apply to surface to ensure complete coverage. For screw cap bottles use the entire contents after opening.

**TECHNICAL DATA:**

Appearance: Clear liquid free from particles. Specific Gravity (20°C): 0.880 to 0.888. Alcohol proportion: 70% +/- 2% v/v.

**STORAGE:**

Store upright in original closed containers away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Material Safety Data Sheet (MSDS).

**SAFE HANDLING:**

Always wear gloves and goggles or face protection. Always read the label and MSDS before use.

**MATERIAL COMPATIBILITY:**

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

**BENEFITS:**

- Low residue.
- Rapid action.
- Bactericidal and Fungicidal
- 3 months validated 'in-use' shelf life
- Contents protected using 'bag-in-bottle' presentation with Premium range.
- Contents protected using filtered trigger presentation with standard range.
- 1L and 5L formats manufactured with WFI quality water.
- Multiple bags for cleanroom Transfer.
- Ideal surface and aseptic transfer Disinfectants.
- Manufactured to GMP.

**TECHNICAL USER INFORMATION:**

InSpec DE is a sterile ready to use disinfectant. This is a 70% v/v alcohol solution in WFI quality water.

InSpec DE is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment. The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (at 25-45kGy) to give a sterility assurance level SAL of 10<sup>-6</sup>.

InSpec DE is supplied with Certificates of Conformity, Analysis, Irradiation Sterility and Endotoxin.

All formats of DE have a stabilised 24 month shelf life.

Microbiological Minimum Efficacy Testing Contact			
EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes

InSpec™ DE Product			
Product	Pack Size	Case Size	Product Code
InSpec™ DE 1L Premium	1L trigger spray	6 x 1L	DEWFI10-1LS
InSpec™ DE 5L Screw Cap	5L screw cap	2 x 5L	DEWFI10-5LS
InSpec™ DE 500ml Premium	500ml trigger spray	8 x 500ml	DEWFI11-500MLS

**Vietnam Official Distributor:**

**Decon Solutions Co., LTD**

No 62 Road 64, Thanh My Loi Ward, District 2, HCM City, Vietnam  
**H:** 09 0890 1955    **F:** +84 28 6287 3181    **E:** [info@deconsolutions.com.vn](mailto:info@deconsolutions.com.vn)  
**T:** +84 28 6287 3180 - 38 143 585    **W:** [www.deconsolutions.com.vn](http://www.deconsolutions.com.vn)