# HA

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Low odour, Safe handling  
 Low residue and Fully effective

**USE INSTRUCTIONS:**

InSpec HA is designed for spraying, wiping and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping.

**TECHNICAL DATA:**

Appearance: Colourless aqueous solution with a characteristic mild chlorine odour. pH 6.5 to 7.5.

**STORAGE:**

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Material Safety Data Sheet (MSDS).

**SAFE HANDLING:**

Always wear gloves and goggles or face protection. Always read the label and MSDS before use.

**MATERIAL COMPATIBILITY:**

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

**BENEFITS:**

- One of the safest and fully effective sporicides. All formats are non hazardous
- Very low residue, corrosion and odour
- Stabilised solution - 12 months shelf life
- 3 months validated 'in-use' shelf life
- Contents protected using 'bag-in-bottle' presentation with premium range.
- 1L and 5L formats available as Ready to Use.
- Multiple bags for cleanroom transfer
- Ideal rotational partner for InSpec QT and AN.
- Manufactured to GMP.

**TECHNICAL USER INFORMATION:**

InSpec HA is a sterile sporicide available ready to use.

InSpec HA is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and aseptically filled into pre-gamma irradiated packaging (25-45kGy) and tested to give a sterility assurance level SAL of  $10^{-6}$ .

InSpec HA is supplied with Certificates of Conformity, Analysis, Irradiation and Sterility.

All formats of HA have a stabilised 12 month shelf life.

For screw cap bottles use the entire contents after opening.

Microbiological Minimum Efficacy Testing Contact			
EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes
EN 13704	1 minute ( Bacillus Subtilis )		

InSpec™ HA Product			
Product	Pack Size	Case Size	Product Code
InSpec™ HAIL Trigger	1L trigger spray	6x 1L	HAPW50-1LS
InSpec™ HA 5L Screw Cap	5L screw cap	2 x 5L	HAPW50-5LS
InSpec™ HA Burstable Pouch	15 wipes/pouch	20 pouches	HAPW50BP15S

**Vietnam Official Distributor:**

**Decon Solutions Co., LTD**

No 62 Road 64, Thanh My Loi Ward, District 2, HCM City, Vietnam  
**H:** 09 0890 1955    **F:** +84 28 6287 3181    **E:** [info@deconsolutions.com.vn](mailto:info@deconsolutions.com.vn)  
**T:** +84 28 6287 3180 - 38 143 585    **W:** [www.deconsolutions.com.vn](http://www.deconsolutions.com.vn)